# FZ / FX / FZ-WP / FX-WP Series

Quick Start Guide  
FZ/FX/FZ-WP/FX-WP Series  
Precision Balances

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A&D Company, Ltd.

## WARNING DEFINITIONS

The warning described in this document has the following meaning:

 <b>CAUTION</b>	A potentially hazardous situation which, if not avoided, may result in personal injury or property damage.
--	--

© 2023 A&D Company, Limited. All rights reserved.

No part of this publication may be reproduced, transmitted, transcribed, or translated into any language in any form by any means without the written permission of A&D Company, Limited.

The contents of this publication and the specifications of the instrument covered by this publication are subject to change for improvement without notice.

Product names and company names mentioned in this document are trademarks or registered trademarks of their respective companies in Japan or other countries and regions.

## Contents

Quick Start Guide – FZ/FX/FZ-WP/FX-WP Series	<b>EN</b>
Precision Balances .....	4
Guide de démarrage rapide – Séries FZ/FX/FZ-WP/FX-WP	<b>FR</b>
Balances de précision .....	15
Kurzanleitung – Baureihe FZ/FX/FZ-WP/FX-WP	<b>DE</b>
Präzisionswaagen .....	27
Guía de inicio rápido – Serie FZ/FX/FZ-WP/FX-WP	<b>ES</b>
Básculas de precisión .....	39
Guida introduttiva rapida – Serie FZ/FX/FZ-WP/FX-WP	<b>IT</b>
Bilance di precisione .....	52
快速入門指南– FZ/FX/FZ-WP/FX-WP 系列	<b>TC</b>
精密天平 .....	64

# 1. Quick Start Guide

Thank you for purchasing an A&D FZ / FX / FZ-WP / FX-WP series precision balance. This quick start guide describes how to install the FZ / FX / FZ-WP / FX-WP series balance, and it covers the balance's basic functions and main operations.

For further information about the FZ / FX / FZ-WP / FX-WP series balance, please refer to the FZ / FX / FZ-WP / FX-WP Series Instruction Manual.

The FZ / FX / FZ-WP / FX-WP Series Instruction Manual is available from the A&D website.

<https://www.aandd.jp>

# 2. Assembly and Installation

An FZ / FX / FZ-WP / FX-WP series balance is a precision instrument and should be carefully unpacked.

It is advisable to store the packing materials so that they can be used when transporting the balance.


The contents of the package vary depending on the model. Refer to the illustration of the packing contents and make sure that everything is included.

## ⚠ CAUTION

- Please confirm that the AC adapter type is correct for your local voltage and receptacle type.
- Use the dedicated AC adapter specified for the balance.
- Do not connect the included AC adapter to other devices.
- If you use the wrong AC adapter, the balance and other equipment may not work properly.


### FZ / FX, 0.0001 g models

A large breeze break is provided as standard for all the models.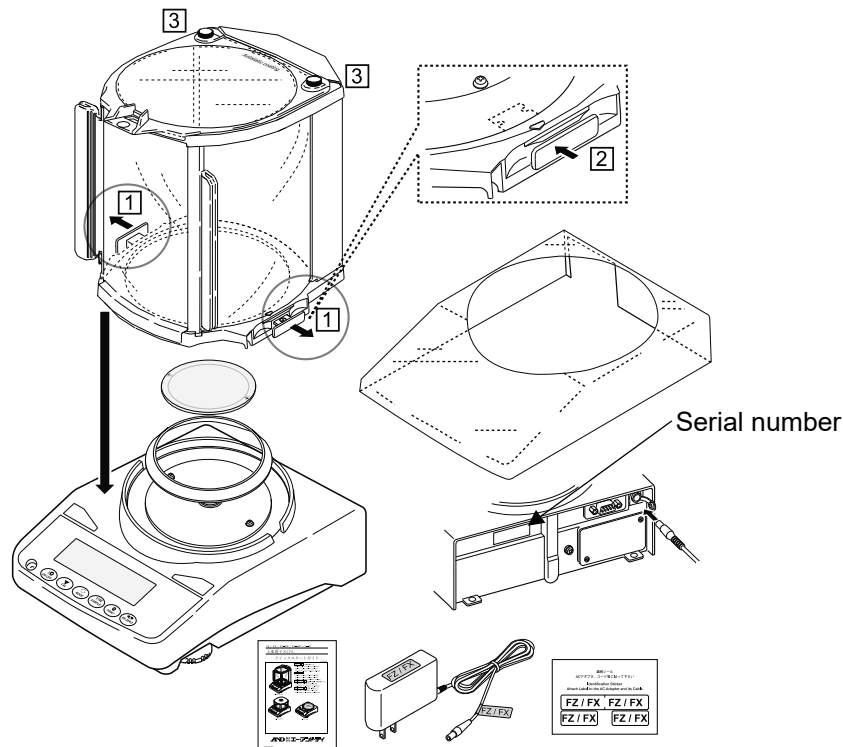


**FZ / FX, 0.001 g models / 0.01 g models**

A small breeze break is provided as standard for all the FZ models and FX 0.001 g models.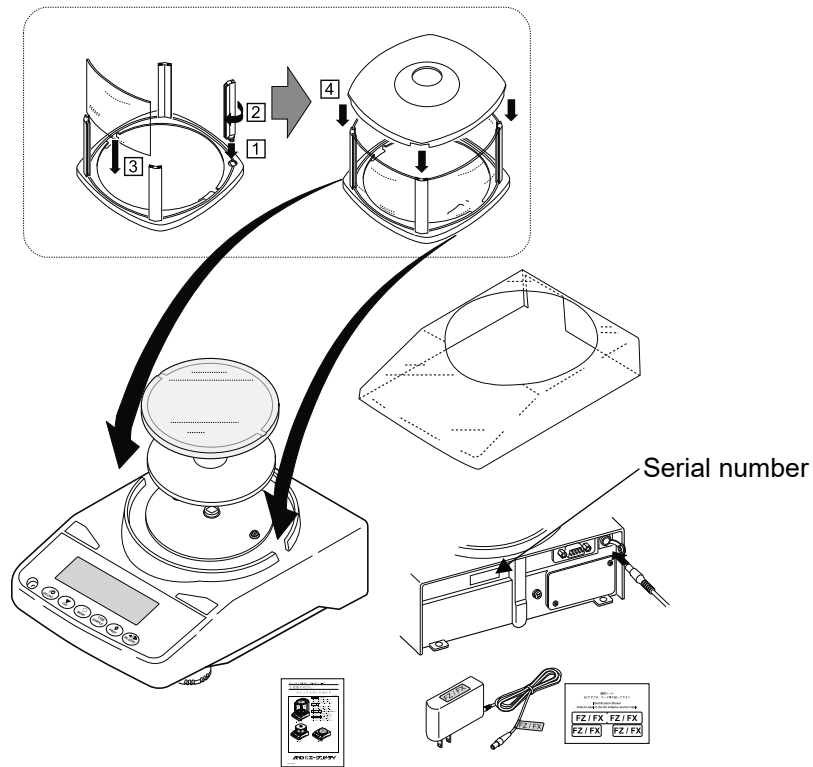

**FZ-WP / FX-WP, 0.001 g models / 0.01 g models**

A small breeze break is provided as standard for all the FZ-WP models and FX-WP 0.001 g models.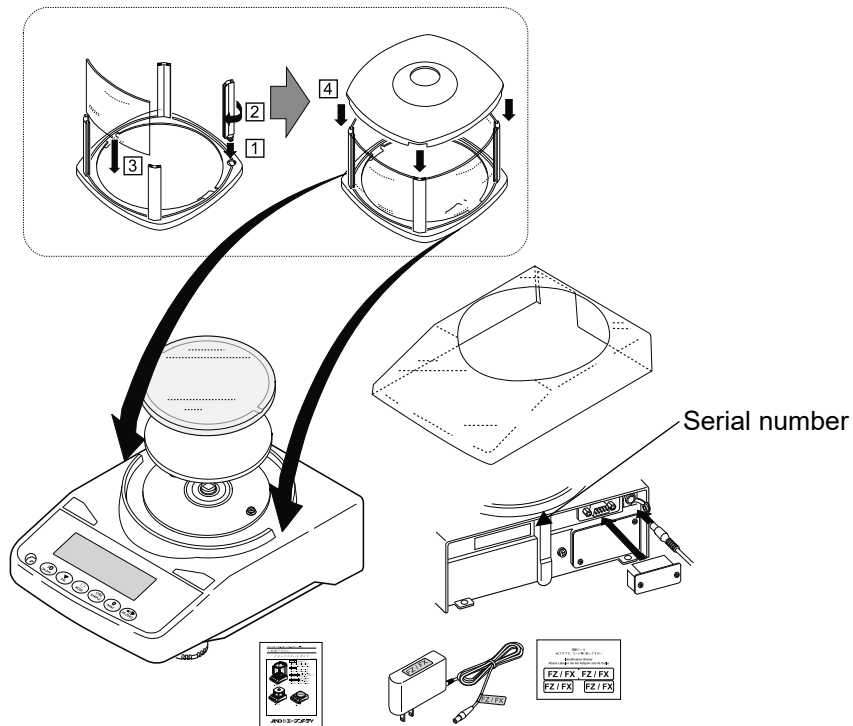


## 2.1. Installation Considerations, Preparation, and Precautions

Prepare the following installation conditions to bring out the full performance of the balance.

- Install the balance in an environment where the temperature and humidity are not excessive. The best operating temperature is about  $20^{\circ}\text{C} \pm 2^{\circ}\text{C}$  at about 45% to 60% RH relative humidity.
- Install the balance where it is free of dust.
- The weighing table should be solid. (An anti-vibration table or stone table is ideal.)
- Place the balance on a horizontal table, and make sure that it is not tilted.
- Install the balance in a stable location, avoiding vibration and shock. The corners of rooms on the first floor are best, as they are less prone to vibration.
- Install the balance where it is not affected by heaters or air conditioners.
- Install the balance where it is not exposed to direct sunlight.
- Install the balance away from equipment that produces magnetic fields.
- Level the balance with the leveling feet and bubble spirit level. Refer to "How to adjust the level of the balance" on the next page.
- Be sure to warm up the balance before use for at least half an hour, or at least an hour for the 0.0001 g model, with the AC adapter connected to the power supply.
- Adjust the sensitivity of the balance before using it for the first time or after having moved it to another location so that accurate weighing can be performed.

For details, refer to "7. Sensitivity Adjustment / Calibration Test" in the separate instruction manual.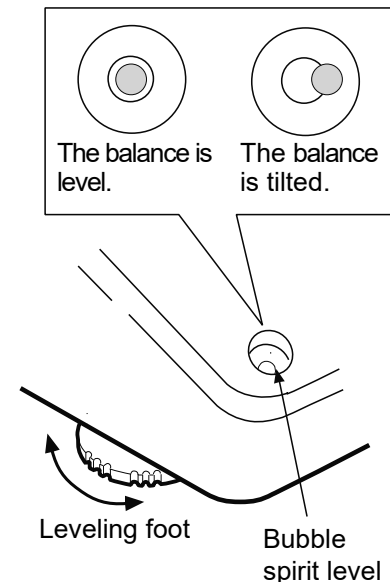


About the FZ-WP / FX-WP series

- The dustproof and waterproof level of the balance is equivalent to IP65, and its second digit, "5", corresponds to "having no harmful influence by receiving direct jet of water". Washing with strong water pressure or submersion in water may cause water to enter the balance and cause a malfunction.
- When installing and using the balance under conditions requiring dustproof and waterproof performance, make sure that the AC adapter plug is fully inserted into the AC adapter jack and that the terminal cover is attached to the RS-232C interface or the waterproof RS-232C cable (AX-KO2737-500) is used.
- If the RS-232C terminal cover is removed or the waterproof RS-232C cable (AX-KO2737-500) is not used, protection against dust and water is not provided.

### CAUTION

Do not install the balance where flammable or corrosive gas is present.

# How to adjust the level of the balance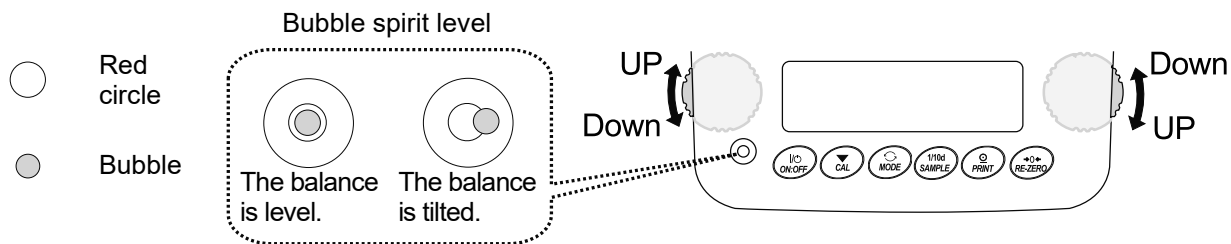

**When the bubble is off to the left.**

Turn the leveling foot on the front right in the clockwise direction.

**When the bubble is off to the right.**

Turn the leveling foot on the front left in the clockwise direction.

**When the bubble is off to the backward position.**


Turn both leveling feet on the front in the clockwise direction at the same time.

**When the bubble is off to the forward position.**


Turn both leveling feet on the front in the counterclockwise direction at the same time.

### 3. Display and Key Panel (Basic Operation)

Lit display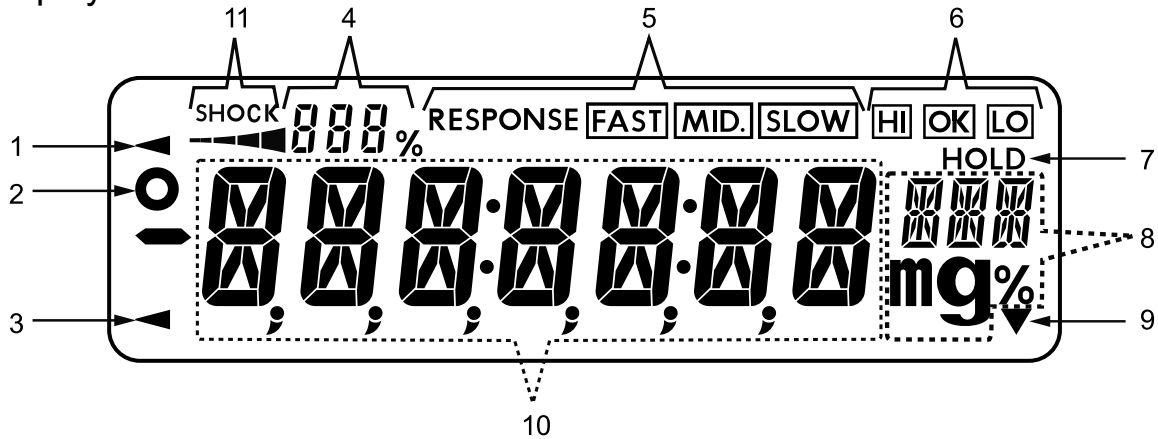

Blinking display

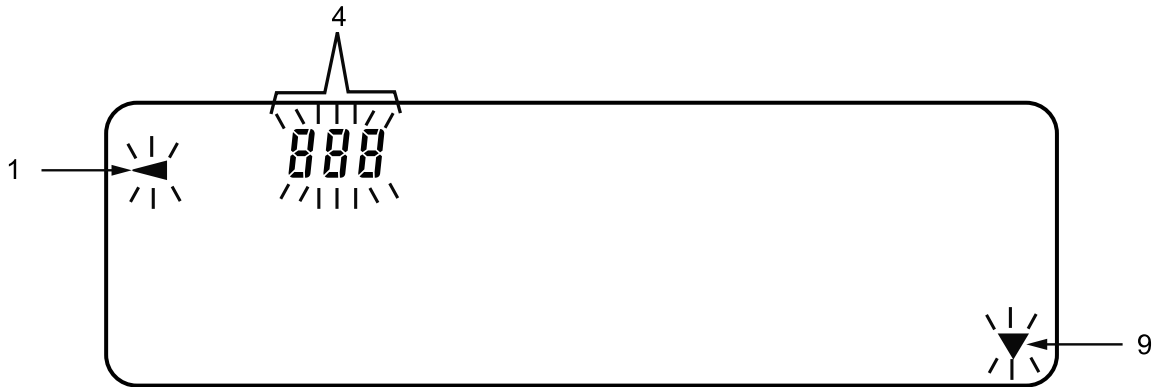No.	Name
1	Processing indicator
2	Stabilization indicator
3	Standby indicator
4	Number of statistical data (Statistical calculation mode) Displays the weight data relative to the weighing capacity, in percentage, in weighing mode (Capacity indicator) Function table setting value display
5	Response indicator (lights for 30 seconds after start of weighing)
6	Comparator indicators
7	Animal weighing indicator
8	Unit display Display for the function table settings
9	Interval output mode indicator (When lit: In standby mode. When blinking: Running)
10	Displays the weighing value or setting item name
11	Shock indicators (ISD)

## Key operation

Key operations affect how the balance functions. Normal key operation during measurement is “press and release the key immediately” or “press the key”. Please do not “press and hold the key (for 2 seconds)” unless required.

 <p>Press the key (Press and release the key immediately.)</p>	 <p>Press and hold the key (for 2 seconds)</p>
---	---

Key	When pressed and released	When pressed and held (for 2 seconds)
	Turns the display on / off. The standby indicator is displayed when the display is turned off. The weighing mode is enabled when the display is turned on. The <b>ON:OFF</b> key is active anytime. Pressing the <b>ON:OFF</b> key during operation will always turn off the display.	
	In the weighing mode, turns the digit for readability on / off. In the counting or percent mode, enters the sample storing mode.	Enters the function table mode.
	Switches the weighing units stored in the function table.	Enters the mode for changing the weighing speed.
	Enters the mode for sensitivity adjustment using the internal weight (FZ / FZ-WP series only). In the function table setting mode, the operation is canceled.	Enters the mode for sensitivity adjustment using an external weight (with your weight).
	When stable, outputs the weighing data according to the function table. In the function table, confirms the operation.	No function at the time of shipment. Outputs "Title block" and "End block" for GLP, GMP, etc. report by changing the function table.
	Sets the displayed value to zero.	

## 4. Sensitivity Adjustment

### 4.1. Sensitivity Adjustment Using the Internal Weight (FZ / FZ-WP Series Only)

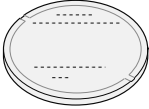

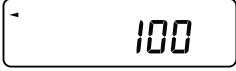
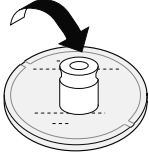
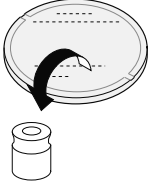
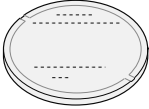
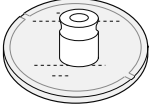
Sensitivity adjustment using the internal weight can be performed with one key press.

Step	Description	Display / Key	Weighing
1	Connect the AC adapter and warm up the balance for at least 30 minutes, or at least an hour for 0.0001 g models, with nothing on the weighing pan.		
2	Press the <b>CAL</b> key. The balance displays	 	 or 
3	The balance automatically performs sensitivity adjustment using the internal weight. Do not apply vibration and the like to the balance.	 	
4	The display automatically returns to weighing mode.	 	



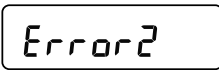
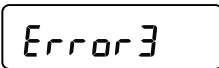
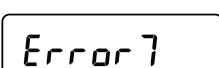
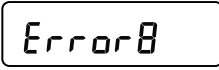
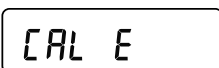
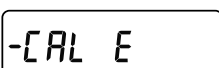
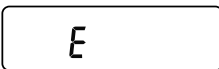
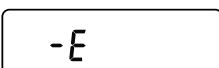
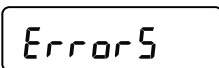
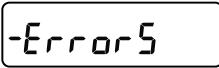
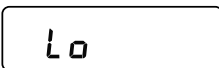
### 4.2. Sensitivity Adjustment Using an External Weight


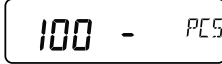
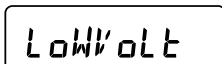
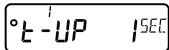
This function performs sensitivity adjustment of the balance using your external weight.

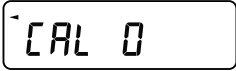Step	Description	Display / Key	Weighing
1	Connect the AC adapter and warm up the balance for at least 30 minutes with nothing on the weighing pan.		
2	Press and hold the <b>CAL</b> key for 2 seconds until  is displayed. Holding down the key switches the items (about every 2 seconds).	 Press and hold (for 2 seconds)   The display switches.	

Step	Description	Display / Key	Weighing
3	Release the key when  is displayed.	  Release the key.	
4	To set the weight value, refer to "7.5. Setting the value of the weight" in the separate instruction manual. For available weights, refer to "7. Sensitivity Adjustment / Calibration Test" in the separate instruction manual. If you do not want to change the weight, proceed to step 5.		
5	Make sure that nothing is on the weighing pan and press the  key to weigh the zero point. Do not apply vibration and the like to the balance.		
6	Place the external weight on the weighing pan and press the  key to weigh the weight. Do not apply vibration and the like to the balance.	  	
7	Remove the external weight from the weighing pan.		
8	The display automatically returns to weighing mode.		
9	Place the external weight again and check that the set value is $\pm 2$ digits. If it is out of range, pay attention to the surrounding environment and start from step 1.		

## 5. Error Codes

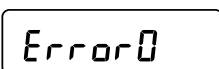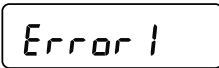Display	Code	Description
		<b>Internal error</b> If this error is displayed continuously, contact your local A&D dealer to repair the balance.
	EC, E11	<b>Stability error</b> Weighing value is unstable and "zero display", "sensitivity adjustment", etc. cannot be executed. Check around the pan. Refer to "2.7. Precautions during use for more accurate weighing" in the separate instruction manual. Improve the environment of the installation location (vibration, drafts, static electricity, etc.). To return to weighing mode, press the  key.
		<b>Out of the setting range</b> The value to be set exceeds the setting range. Enter again within the setting range.
		<b>Malfunction of the internal memory element of the balance</b> If this error continues to be displayed, please request a repair.
	EC, E17	<b>Internal weight error (FZ / FZ-WP series only)</b> The internal weight application mechanism does not function properly. Perform the operation from the beginning. If the error does not improve even after performing the operation again, please request a repair.
		<b>Abnormality in the internal memory data of the balance</b> If this error continues to be displayed, please request a repair.
	EC, E20	<b>Sensitivity adjustment weight error (Positive value)</b> The sensitivity adjustment weight is too heavy. Check around the pan. Check the mass value of the weight. To return to weighing mode, press the  key.
	EC, E21	<b>Sensitivity adjustment weight error (Negative value)</b> The sensitivity adjustment weight is too light. Check around the pan. Check the mass value of the weight. To return to weighing mode, press the  key.
		<b>Overload error</b> A sample beyond the balance weighing capacity has been placed on the pan. Remove the object from the pan.
		<b>Weighing pan error</b> The weighing value is too light. Check that the weighing pan is installed correctly. Set the weighing pan correctly. Perform sensitivity adjustment.
		<b>Overload error</b> A sample beyond the balance weighing capacity has been placed on the pan. Remove the object from the pan. If the error cannot be resolved, contact your local A&D dealer.
		<b>Weighing pan error</b> The weighing value is too light. Check that the weighing pan is installed correctly. Set the weighing pan correctly. If the error cannot be resolved, contact your local A&D dealer.
		<b>Sample mass error</b> The sample is too light to be stored as a sample mass for the counting mode or percent mode. The sample cannot be used.

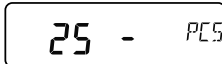Display	Code	Description
  		<b>Unit weight error</b> The sample mass for the counting mode is too light. Storing and using it for counting may cause a counting error. Add samples to reach the specified number and press the <b>PRINT</b> key. Although pressing the <b>PRINT</b> key without adding samples will put the balance in counting mode, make sure to add samples for accurate counting.
		<b>Clock battery error</b> The clock backup battery has been depleted. Press any key and set the time and date. Even if the clock backup battery is depleted, the clock and calendar function works normally as long as the balance is powered with the AC adapter. If this error appears frequently, contact your local A&D dealer for repair.
		<b>Power supply voltage fault</b> The voltage supplied from the AC adapter is abnormal. Check that the AC adapter is the one supplied with the balance.
	EC, E00	<b>Communications error</b> A protocol error occurred in communications. Check the format, baud rate, etc.
	EC, E01	<b>Undefined command error</b> An undefined command was found. Check the transmitted command.
	EC, E02	<b>Not ready</b> The received command cannot be executed: (e.g.) Q command was received when not in weighing mode. (e.g.) Q command was received while re-zeroing. Adjust the delay time to transmit a command.
	EC, E03	<b>Timeout error</b> With the timeout parameter set to  , there was a wait time of approximately 1 second or more while receiving command characters. Check the communication.
	EC, E04	<b>Character length error</b> The number of characters in the received command has exceeded the limit. Check the command to transmit.
	EC, E06	<b>Format error</b> The description of the received command is incorrect: (e.g.) The number of digits of numerical values is incorrect. (e.g.) There are alphabet characters among the numerical values. Check the transmitted command.
	EC, E07	<b>Parameter setting error</b> The value of the received command has exceeded the allowed value. Check the setting range of the numerical value of the command.
Other error code		If an error other than the above is displayed, or if any of the above errors cannot be resolved, please contact your local A&D dealer.

## 6. Specifications

	FZ / FX Series			
	104	154	254	254D
Weighing capacity	102 g	152 g	252 g	
Maximum display	102.0084 g	152.0084 g	252.0084 g	252.008 g
				62.0009 g *1
Readability	0.0001 g			0.001 g
				0.0001 g
Repeatability (Standard deviation)	0.0001 g		0.0002 g (200 to 250 g) 0.0001 g (0 to 200 g)	0.0005 g
				0.0001 g
Linearity	±0.0002 g		±0.0003 g	±0.001 g
				±0.0003 g

\*1 Precision range weighing is possible even with a heavy tare. (Smart Range Function)

	FZ/ FX Series FZ-WP / FX-WP Series *3							
	123 123WP	223 223WP	323 323WP	523	1202 1202WP	2202 2202WP	3202 3202WP	5202
Weighing capacity	122 g	220 g	320 g	520 g	1220 g	2200 g	3200 g	5200 g
Maximum display	122.084 g	220.084 g	320.084 g	520.084 g	1220.84 g	2200.84 g	3200.84 g	5200.84 g
Readability	0.001 g				0.01 g			
Repeatability (Standard deviation)	0.001 g				0.01 g			
Linearity	±0.002 g				±0.02 g			
Accuracy after sensitivity adjustment using the internal weight *2 (FZ / FZ-WP series only)	±0.010 g				±0.10 g		±0.15 g	

\*2  Accuracy after sensitivity adjustment using the internal weight, with the accessory small breeze break installed and in good ambient conditions within the temperature range of 10°C to 30°C with no abrupt changes in temperature or humidity, no drafts, no vibration and no effect by magnetic fields or static electricity.

With the FZ / FZ-WP series balance, the environment and aging may cause internal weight value errors. The balance requires regular maintenance and adjustment using an external weight.

\*3 The FZ-WP / FX-WP series balances have dustproof and waterproof performances (IP65 compliant).

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### **A&D AUSTRALASIA PTY LTD**
### **A&D KOREA Limited**
### **A&D Instruments India Private Limited**