


Instruction Manual for AD-1676 Tabletop Breeze Break (M)

1. Features


- The tabletop breeze break protects the balance from breezes due to air conditioning and body motion, thereby reducing balance weighing error.
- This size of tabletop breeze break is most suitable for use with the BM series.
- The transparent panel assembly consists of antistatic plastic material, which protects the balance from static electricity.
- The hole on the back panel can be used to pass the RS-232C and AC adapter cables through.

2. Specifications

- Weight : Approx. 4 kg
- Size : 356 (Width) x 514 (Depth) x 350 (Height) mm
- Material : Transparent panel: Antistatic plastic
Rear panel / Frame: Aluminum
- Applicable models (BM series balance)
- * We recommend installing the tabletop breeze break (M) in a position where the display of the balance can be seen easily and the balance can be operated properly.
- * To minimize the effect of ambient drafts on measurement, open the side door A only as far as necessary.

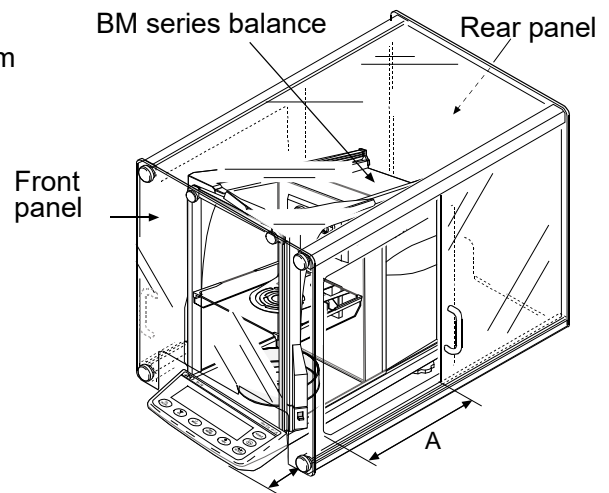

3. External dimensions

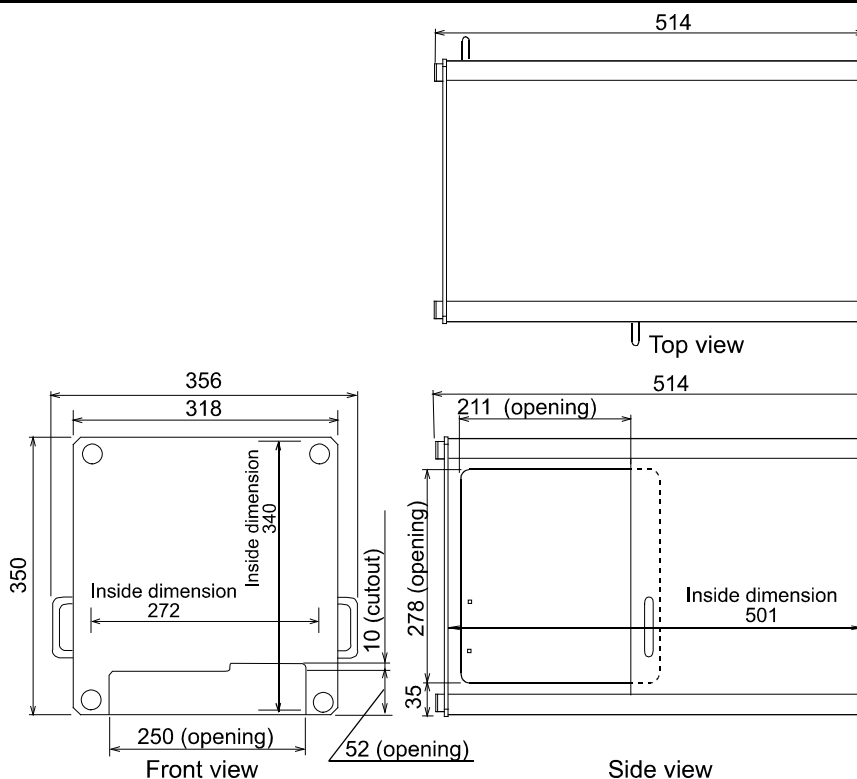

Caution

- Do not use organic solvents to clean the tabletop breeze break as that may damage the transparency and antistatic properties of the panels.
- Use a soft lint free cloth dampened with warm water and a mild detergent when cleaning the tabletop breeze break.