
HJ-150 Compact Scale

Operating Instructions

Cautions

- ❑ Avoid installing / leaving the scale in direct sunlight / high humidity.
- ❑ Place the scale on a flat and solid surface in a place free from wind.
- ❑ Avoid vibration, strong electric or magnetic fields and rapid changes in temperature or humidity.
- ❑ Avoid using the scale near other electronic equipment.
- ❑ Do not mix the battery types. Replace all of the batteries at same time.
- ❑ If the scale is not to be used for a long period of time, remove the batteries.
- ❑ Avoid over / shock load that could cause damage to the scale.
- ❑ Keep the protective cover attached when you carry the scale.
- ❑ Wait for more than 6 seconds to let the scale warm up after turning on the power.
- ❑ When the location to be used or the ambient environment has changed, calibrate the scale before use.
- ❑ When the display shows $L \square$, or continues to show $\square\square\square\square$ more than 30 seconds after turnig on, replace the batteries with new one.

Specifications

Capacity: 150 x 0.1 g / 5.29 x 0.01 oz

Linearity: ± 0.2 g Repeatability: 0.1 g (Standard deviation)


Operating environment: 5°C ~ 40°C (41°F ~ 104°F), less than 85% RH

Display: 7 segment LCD with backlight (Character height 8 mm)

Power: 2 x "AAA" size batteries, approx. 80 hours with alkaline batteries.

Pan size: 60 (W)x 64 (D) mm / 2.36 (W) x 2.52 (D) in.

Accessories: Instruction manual, 100g weight, 2 monitor batteries

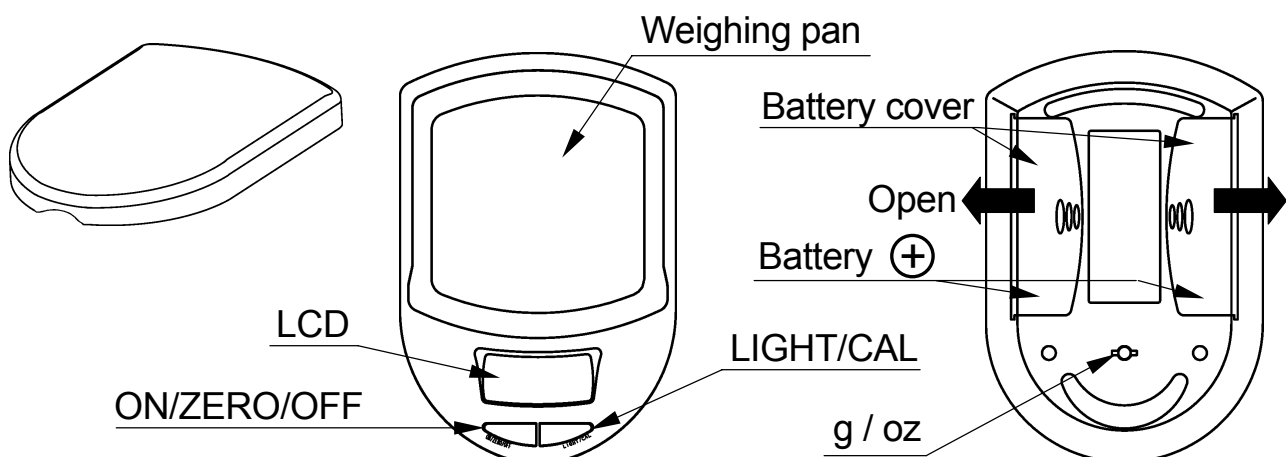

Switches

ON/ZERO/OFF Switch:

Press to turn the scale on. In the operation mode, press briefly to set the display to zero. Press and hold approximately 3 seconds to turn the scale off.

LIGHT/CAL Switch:

In the operation mode, press briefly to turn the LCD backlight on. Press and hold approximately 3 seconds to enter the calibration mode.

g / oz Switch:

In the operation mode, press the switch located on the bottom of the scale to change the weighing unit between “g” and “oz”.

Installing / Replacing batteries

Remove the battery covers and insert 2 AAA size batteries into each housing, taking care with the polarities. Attach the battery covers again.

- Keep the protective cover attached when installing and/or replacing batteries, to prevent excessive force from being added to the weighing pan.

Weighing procedures

1. Turn on / off the scale

Press the **ON/ZERO/OFF** key to turn on the scale. The scale shows all segments for a few seconds and displays zero. The backlight will turn on for about one minute. Press and hold approximately 3 seconds to turn the scale off.

- When powered on and unstable weighing continues more than 15 seconds, **8.8.8.8** display goes to **E** and the scale turns off.


2. Select the weighing unit

Press the **g / oz** switch to select a weighing unit from “g” and “oz”.

3. Weighing

Verify the reading is zero. If not, press the **ON/ZERO/OFF** key to zero the display. If you use a container for weighing, place an empty container on the weighing pan and press the **ON/ZERO/OFF** key to zero the display.


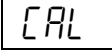
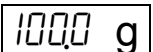
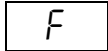
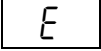
Place the objects to be weighed on the pan or in the container. Wait for the display to become stable and read the weight.

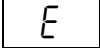 means overload. Remove the objects from the pan immediately.

Automatic power-off function


The automatic power-off function turns power off after approximately 2 minutes of stable display during the object(s) applied. When there is no object applied, it will turn off after 30 seconds.

Calibration

1. Make sure that there is nothing on the weighing pan, press and hold the  key approximately 3 seconds. The display shows . Release the key.
2. The scale memorizes the zero point and automatically displays .
3. Place the weight of 100g on the center of the pan within 10 seconds.
4. After approximately 8 seconds, the scale will display  for success and  for failure. Then, the scale automatically turns off.
5. Turn the scale on again to start the operation.

If calibration fails, check if anything is touching the weighing pan, there is vibration or wind, and the calibration weight is correct.

Dimension

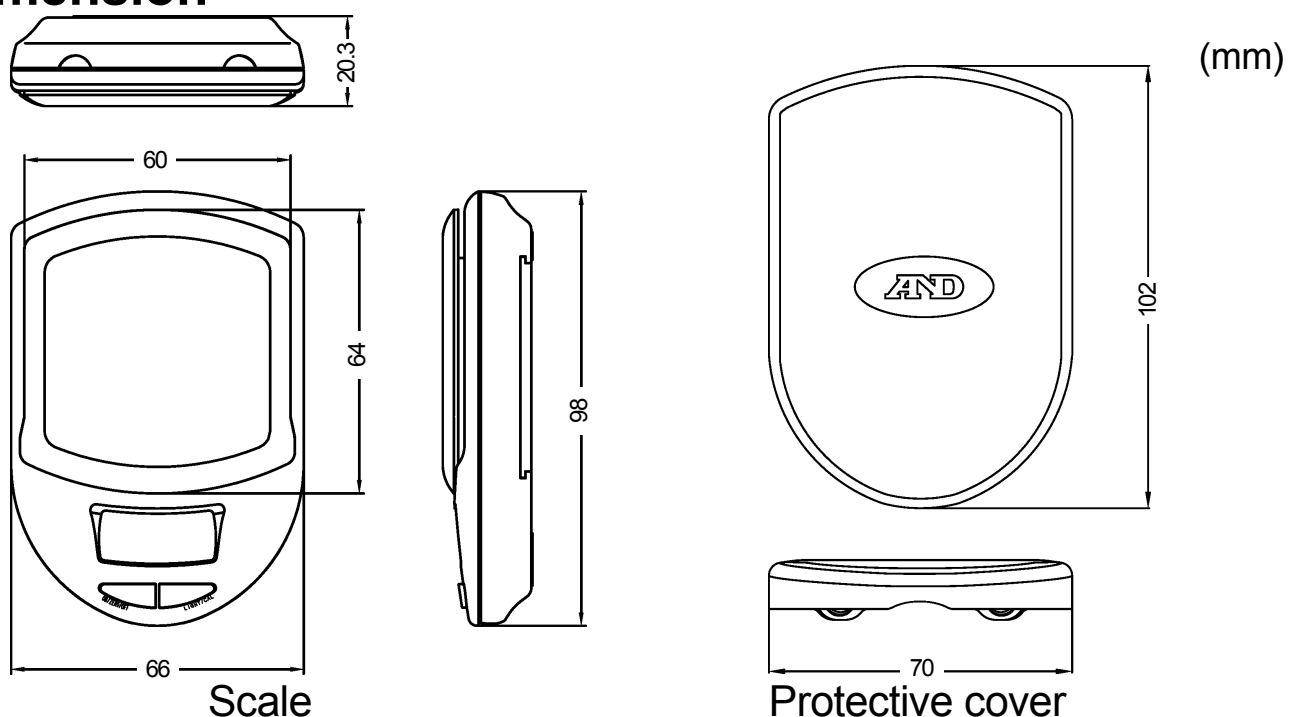



A&D Company, Limited

3-23-14 Higashi-Ikebukuro, Toshima-ku, Tokyo 170-0013 JAPAN
Telephone: [81] (3) 5391-6132 Fax: [81] (3) 5391-6148

A&D ENGINEERING, INC.

1756 Automation Parkway, San Jose, California 95131 U.S.A.
Telephone: [1] (408) 263-5333 Fax: [1] (408) 263-0119

A&D INSTRUMENTS LIMITED <UK Office>

Unit 24/26 Blacklands Way, Abingdon Business Park, Abingdon, Oxfordshire OX14 1DY
United Kingdom
Telephone: [44] (1235) 550420 Fax: [44] (1235) 550485

A&D INSTRUMENTS LIMITED <German Sales Office>

Große Straße 13 b 22926 Ahrensburg Deutschland
Telephone: [49] (0) 4102 459230 Fax: [49] (0) 4102 459231

A&D MERCURY PTY. LTD.

32 Dew Street, Thebarton, South Australia 5031 AUSTRALIA
Telephone: [61] (8) 8301-8100 Fax: [61] (8) 8352-7409

A&D KOREA Limited

8th Floor, Manhattan Bldg. 36-2 Yoido-dong, Youngdeungpo-ku, Seoul, KOREA
Telephone: [82] (2) 780-4101 Fax: [82] (2) 782-4280

A&D RUS CO., LTD.

112 Kuntsevo Block, Vereyskaya st., Moscow, 121357 RUSSIAN FEDERATION
Telephone: [7] (495) 937-33-44 Fax: [7] (495) 937-55-66

A&D Instruments India Private Limited

509, Udyog Vihar, Phase-V, Gurgaon - 122 016, Haryana, India
Telephone: 91-124-4715555 Fax: 91-124-4715599