



HT-CL Series Supplementary Instructions

The following information was changed/added to the HT-CL series instruction manual.

This addition to the instruction manual supports software version P-101 and later.

[Regarding Negative comparison for take away [PP1]]

Take-away check weighing (negative comparison) is the way to compare the negative weight while taking away objects from a container. Set the function [PP1].

Note

To start the take-away check weighing, be sure to use the RE-ZERO key to tare the weight of the container filled with objects. The RE-ZERO key may zero the display, and the scale goes below the zero point by taking out the objects.

Function setting list

Item / Display	Parameter	Description
Normal / Negative comparison [PP]	◆0	Normal comparison
	1	Negative comparison for take-away check weighing

◆ : Initial factory setting

