

Models that can be connected to Mac via USB (WinCT is not supported)

2018/2/6

Direct input to application (Quick USB connection)			
Model	Interface	Remarks	OS Verified
BM	Standard USB interface	No setting required	mac OS X 10.10 Yosemite macOS 10.12 Sierra
GH/HR-i	GH-02		
HR-AZ/A	FXi-02		
MC-1000/6100	GX-02	Internally set to Quick USB NU	
GX-A/GF-A	Standard USB interface		
GX/GF	GX-02	No setting required	
FZ-i/FX-i	FXi-02		
EJ	EJ-02		
SC/SE	SCE-02		

Capture logging data			
Model	Output data	Remarks	OS Verified
Weighing Environment Logger AD-1687	Date + Time, Temperature, Humidity, Barometric Pressure, Vibration and Weighing Data	Read in CSV file *It does not work on Real-Time Transmission Mode	OS X 10.10 Yosemite macOS 10.12 Sierra
Weighing Data Logger AD-1688	Date + Time and Weighing Data	Read in CSV file	

Caution

***Only the weighing values can be received by Mac.**

When data other than weighing value is output from Quick USB, **alphanumeric characters are corrupted.**

The GX-A/GF-A should be set to Quick USB NU mode.

***Quick USB Adapter (AD-8527) does not work on Mac.**