
HL-*i* series Digital Compact Scale

Calibration Instructions

- Warm up the scale for more than 5 minutes. Refer to the instruction manual for how to disable the automatic power-off function, if necessary.

1. Calibration weight

Calibrate the scale with calibration weights equivalent to capacity.

HL-200*i* : 200g ± 0.02g

HL-2000*i*: 2000g ± 0.2g

2. Calibration procedures

- (1) Start with the scale turned off. Press the **ON:OFF** key while pressing both the **RE-ZERO** key and **UNITS** key. **Cal** will be displayed.
 - (2) Press the **RE-ZERO** key to calibrate zero. **Cal 0** will be displayed. Make sure there is nothing on the weighing pan. Wait for the stable indicator to be displayed and press the **RE-ZERO** key. **Cal f** will be displayed in a few seconds.
 - (3) Place the calibration weight on the center of the weighing pan. Wait for the stable indicator to be displayed and press the **RE-ZERO** key. The display will show **end**.
 - (4) Remove the calibration weight and the scale will automatically return to the weighing mode.
 - (5) Place the calibration weight on the center of the weighing pan again and check the calibration has been done correctly.
- If you press the **UNITS** key at step (2), a 4 digit number will be displayed. Press the **UNITS** key again to return to **Cal**.

End of calibration instructions.